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WS 1.6 mm iØ 1.6 – 11.1 mm  
WS 2.4 mm iØ 4.8 + 6.4 mm, (WS = Wandstärke)

## Technical specifications


**Principle:** 1-channel tubing pump, with ball bearings  
3 stainless steel rollers  
adjustable for different tubing types

**Flow-rate:** 0.41– 3600 ml/min  
with 1.6 mm tubing wall thickness  
3.4 – 1500 ml/min  
with 2.4 mm tubing wall thickness

**Pressure:** max. 2.5 bar (36 psi) with tubing up to  
8 mm ID, 1.5 bar (22psi) with larger tubing

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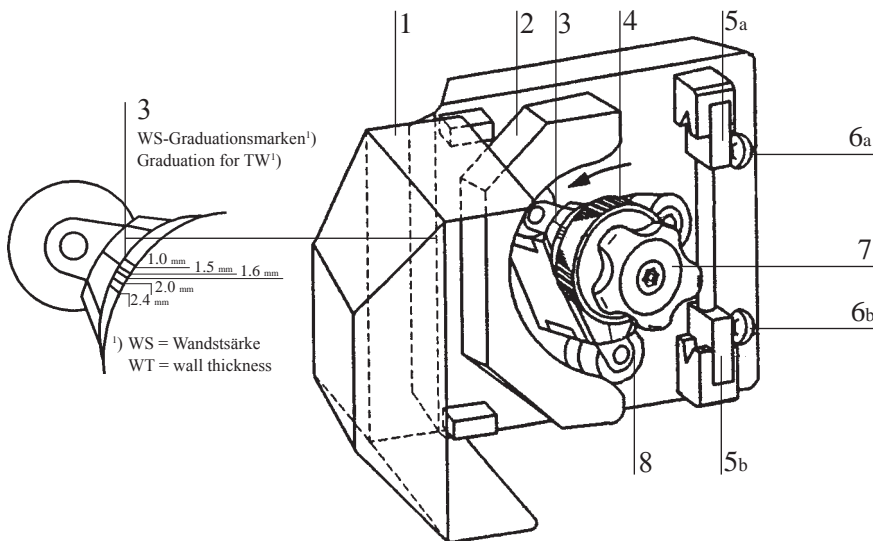## Adjusting the roller pressure

The knurled ring (4) with the 5 graduation marks allows you to adjust the pressure of the rollers onto the tubing as follows:

- Open pumphead cover (1)
- Loosen the star grip (7)  
→ turn in the counter-clockwise direction
- Turn the knurled ring (4) so that the correct graduation mark (depends on the wall thickness, see figure below) is aligned with the notch in the triangular roller-head (3)
- By simultaneous fastening the star grip (7) you lock the pump-rollers in the position required

The particular advantage of this roller adjustment is that the contact pressure of the 3 rollers occurs concentrically from the roller-head shaft to the contour of the tube-bed. This ensures a regular and constant pressure, independently of the position of the roller-head and the wall thickness of the tubing used. This provides a favorable effect on the pump characteristics and the service-life of the tubing.

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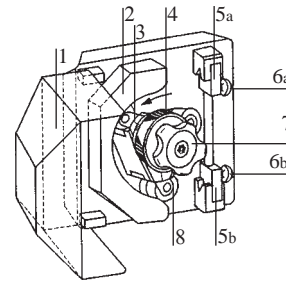
Open the pumphead cover (1)

Load the pump tubing as follows:

- Press the ball handle (6a) down with your thumb
- Insert the tubing with your free hand into the opening of the tubing retainer prism (5a)
- Release the ball handle (6a)
  - ➔ the tubing is now clamped in the tubing retainer prism (5a)
- Insert the tubing into the tube-bed (2) by turning the rotor with the pump rollers (8) in the counter-clockwise direction
  - ➔ the tubing must be well aligned with the tube-bed (2) and centred by the rollers (8)
- Now lift the ball handle (6b) with your thumb and clamp the tubing in the lower tubing retainer prism (5b)

## Start-up procedure

Before switching the drive on, you must close the pumphead cover (1)! Danger of injury!

**Caution:** rotating pumphead creates a pinch and crush hazard.

Before starting the operation of your pump system ensure that the tubing lies correctly in the tube-bed (not twisted) and that the tubing tension is correct. The appropriate tubing tension can be adjusted as follows:

- When pumping in the clockwise direction (\*) press down the upper ball handle (6a)
  - With your other hand pull slightly the tubing so that a weak tension is achieved
  - Release the ball handle (6a)
- 1) When pumping in the counter-clockwise direction, lift the lower ball handle (6b) and adjust the tubing tension.

## MCP pumphead identification

**ID Code 380A.** Must be entered for each of the 4 programs using this pumphead.

Please refer to the drive operating manual.