## Synchronmotor

## Synchronous motor

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Pumpenmodelle

Pump models

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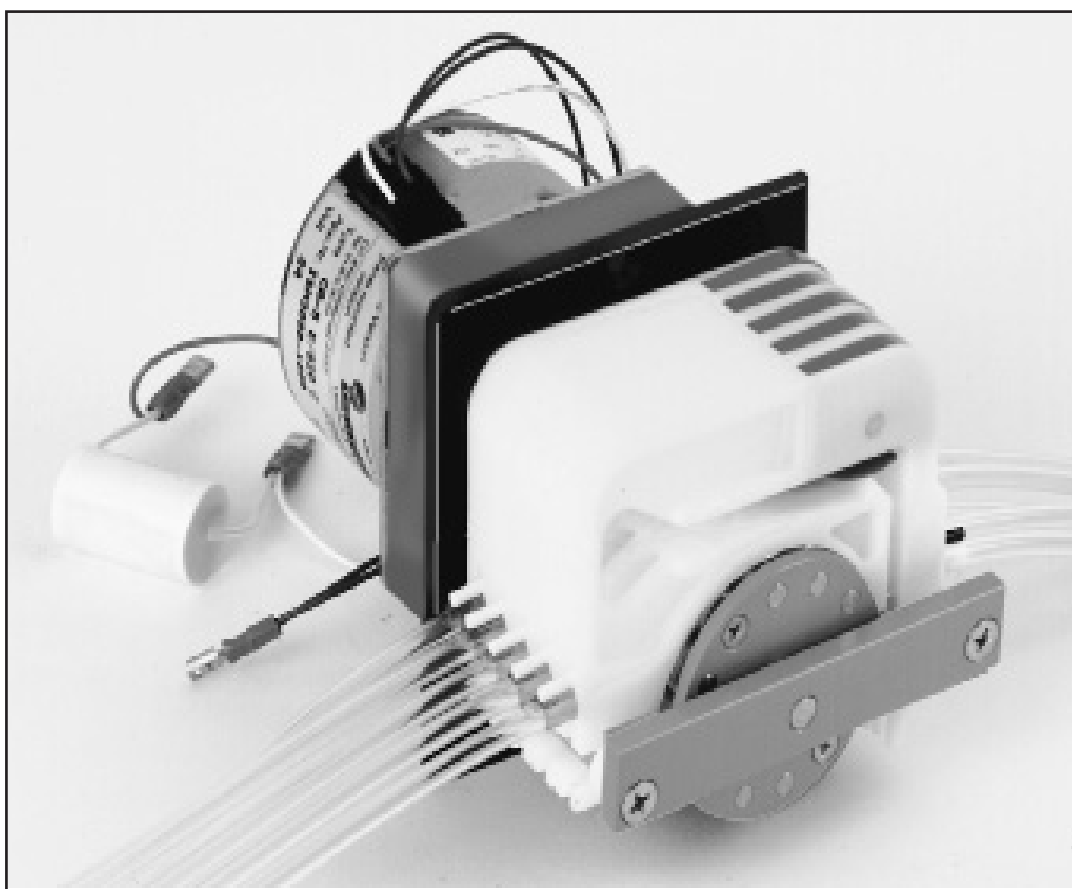
**MS/CA1-E\_**

**MS/CA2-E\_**

**MS/CA3-E\_**

**MS/CA4-E\_**

**MS/CA5-E\_**

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### Contents

- Warranty
- Starting the pump
- Cassettes
- Flow-rates
- Spare-parts list

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## Operating manual for panel mounted tubing pumps

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#### ⚠ Please note

We recommend you read this operating manual carefully.

When operating a pump, certain hazards cannot be excluded.

**ISMATEC® does not take liability for any damage resulting from the use of an ISMATEC® pump.**

ISMATEC® does not admit responsibility for the handling of chemicals.

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#### Safety precautions

ISMATEC® tubing pumps are designed for pumping applications in laboratories and industry. As such it is assumed that Good Laboratory Practice (GLP) and our following recommendations will be observed.

**Caution:** rotating pump-head creates a pinch and crush hazard.

- The circuit between mains supply and pump has to be earthed.
- The pump must not be operated outside the destined operating and environmental conditions.
- The pump **must not** be used:
  - for medical applications on human beings
  - in explosion proof chambers or in the presence of flammable gases or fumes
- The pump must be switched off when cassettes or tubing are inserted or changed.
- Pump-heads consist of revolving parts (rotor of the pump-head). Therefore, the pump must not be operated before the cassettes are fully snapped-in or the tube-bed is completely shut.
- Do not manipulate the pump-head before the pump is switched off and disconnected from the mains.
- Be particularly cautious that no parts of your body such as fingers, long hair, etc. or jewellery, or loose objects such as cables or tubing, etc. can be trapped by the revolving rotor.
- Do not open or remove the housing while the pump is operating.
- Repairs may only be carried out by a skilled person who is aware of the hazard involved.
- For service and repairs carried out by the customer or by third-party companies ISMATEC® denies any responsibility.
- **CE conformity**  
Our standard pump models comply with the applicable EC directives. Within the EC, however, it is not permissible to operate our panel-mounted pumps before the complete system or instrument complies with all the corresponding EC directives.

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#### Warranty

**For all ISMATEC® panel-mounted tubing pumps excluding consumables:**

**from date of delivery: 1 year**

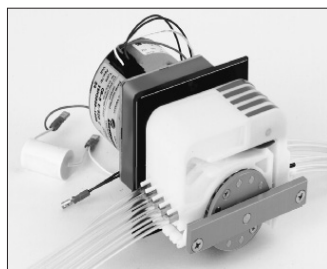
We warrant the perfect functioning of our products, provided they have been installed and operated correctly according to our operating instructions.

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#### Warranty Terms

If production or material faults can be proved, the defective parts will be repaired or replaced free of charge at our discretion. The duration of the warranty is not affected by making a claim for warranty service. Further claims are excluded. Shipping costs for instruments under warranty are charged to the customer.

**Our warranty becomes invalid in case of:**

- improper operation by the user, or if the pump is diverted from its proper use,
- unauthorized modification or misuse by the user or by a third-party,
- improper site preparation and maintenance,
- operation outside environmental and electrical specifications for the product,
- use of third-party software, hardware, interfacing or consumables purchased by the user and which do not comply with our specifications,

#### Product

- Panel-mounted tubing pump MS/CA\_-E\_ with cassettes (pump model as ordered)

Please check the package and its contents for transport damage. If you find any signs of damage, please contact your local ISMATEC® representative immediately.

Complaints can only be accepted within 8 days from receipt of the goods.

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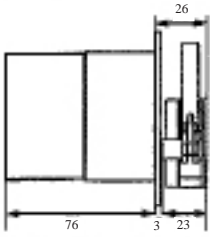
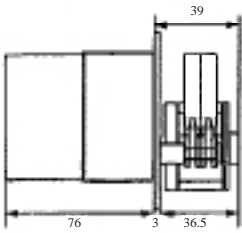
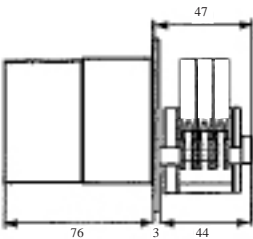
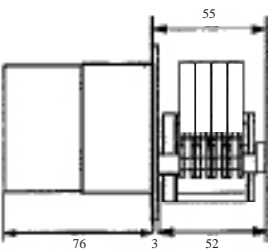
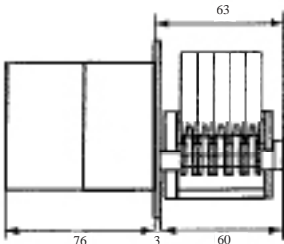
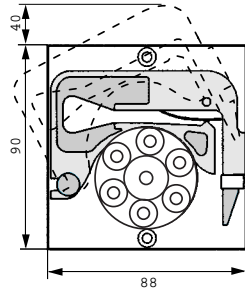
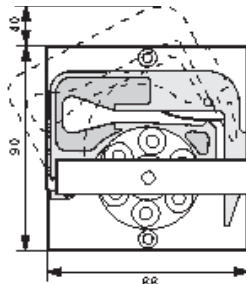
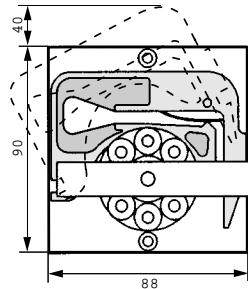
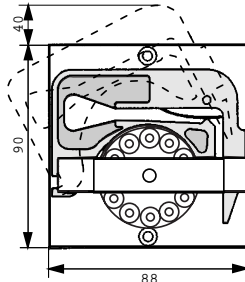
min<sup>-1</sup> x ml pro Umdrehung = ml/min z.B. 20 x 0.43 = 8.6

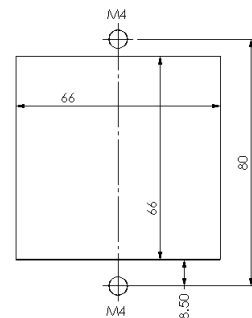
**Example for calculating the flow rate**

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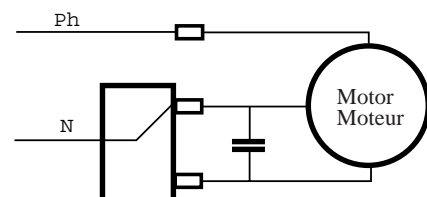
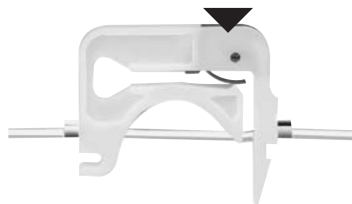
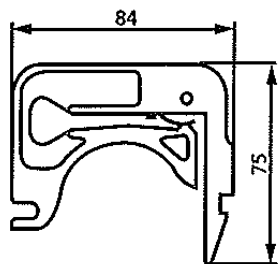
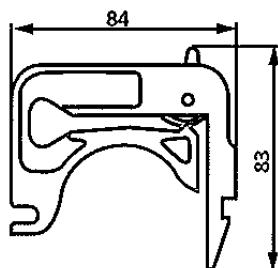
**MS/CA Anpresshebel**  
**MS/CA pressure lever**  
**MS/CA levier de pression**

**3 Color Code Stopper Schläuche**  
**3-stop color-coded tubing**  
**Tubes color-codé à 3 arrêts**

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#### Type of cassettes

**MS/CA Click'n'go**  
Spring-loaded cassette with stainless steel pressure spring providing defined and reproducible occlusion conditions. Independent of tubing diameter and material, the occlusion is automatically correct.

**MS/CA Pressure lever**  
Cassette with pressure lever for adjustable occlusion (may be an advantage for elevated differential pressure conditions).

#### Cassette material

**POM-C**  
(Polyoxymethylene-Copolymer) (Polyacetal, Polyformaldehyde)  
Good chemical resistance to many organic solvents and strong alkaline chemicals. Is affected by strong acids and oxidizing substances.  
UV-stabilized and stable at (dry) temperatures up to 80°C (in continuous use) or 136°C (for a short time).

**PVDF <sup>1)</sup>**  
(Polyvinylidene fluoride)  
Very good chemical resistance to acids and most aliphatic, aromatic and chlorinated solvents.  
Not suitable for long contact with esters, ketones, amines and strong alkaline chemicals.  
Stable to UV and temperatures up to 110°C (continuous use) or 142°C (for a short time)

<sup>1)</sup> only cassettes with pressure lever

#### Pump tubing

These cassette pumps take tubing with 3 color-coded collars. Thanks to these collars the tubes are always inserted under the same conditions, which results in a high repeatability of the flow rate. What is more, the different color codes help to easily recognize the various tubing diameters.

Tubing with 3 collars can be used twice. Thanks to this advantage the service life of the tubes can be doubled.

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#### Type of cassettes

- 1 Remove the cassette by slightly pressing the fixing-tongue and lifting it simultaneously.
- 2 - **Click'n'go Cassettes**  
Insert the tubing (with 3 color-coded collars).  
- **Cassettes with pressure lever**  
Push the pressure level down (it must be relaxed). Insert the tubing (with 3 color-coded collars).
- 3 - **Click'n'go Cassettes**  
Reinsert the cassette(s) into the roller-head.  
- **Cassettes with pressure lever**  
Reinsert the cassette(s) into the roller-head. Lift the pressure lever step by step until the medium is sucked in and a continuous flow is achieved (no further step than necessary!)

#### Please note

- ➔ The cassettes allow the user to insert tubes with different diameters and materials on the same roller-head.
- ➔ When the pump is idle, release all cassettes at the fixing-tongue (2).

⚠ When using new tubing for the first time, it may occur that, depending on the tubing used (hardness and diameter), the pump cannot be primed and, hence, does not deliver the liquid. If that is the case we recommend you to wet the tubing and to run the pump with the tubing inserted for about 15 to 30 minutes.

- ➔ For applications with a high differential pressure (>1 bar), we recommend the use of cassettes with pressure lever.

#### Chemically resistant tubing

**Tygon® MHLL**  
Chemically resistant to acetone methylethylketone and other highly aggressive solvents.

Please also refer to our detailed tubing documentation.

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