REGLO-E Digital, REGLO-E-CPF Digital
Installation und elektrische Anschlüsse
Installation and electrical connections / Installation et connexions électriques

REGLO-E Digital
Einbaupumpen

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Important information

■ **Safety**

The installation of this panel-mounted pump must be carried out by a skilled person who is aware of the potential hazard involved in handling electrical components.

Caution: rotating pumphead creates a pinch and crush hazard.

■ **Wiring**

See figure on front page as well as the attached wiring diagram.

■ **Active cooling system**

Depending on the installation and the operating time an active cooling system (ventilator) must be installed.

■ **Connecting cable**

Shielded cables are to be used between control unit and drive. The shield must be connected to ground on both ends. The maximum length of the cable checked and approved by ISMATEC® is 2 metres. Depending on the situation, longer cables can be used which, however, may cause malfunction. For more information, please do not hesitate to contact us.

■ **CE conformity**

Our standard pump models comply with the applicable EC directives.

Within the EC, however, it is not permissible to operate our panel-mounted pumps before the complete system or instrument complies with all the corresponding EC directives.

Technical documentation

- Please refer to the enclosed
- dimensioned diagrams
 - wiring diagrams
 - operating instructions for:
 - fuse ratings and voltage
 - DIP-switch settings
 - how to operate the pump

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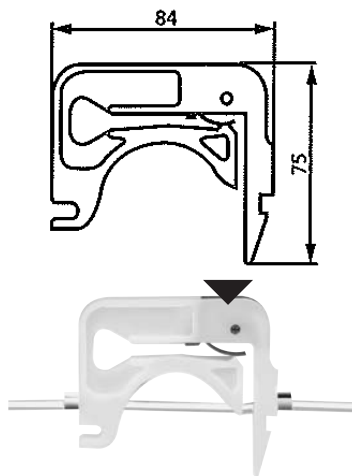
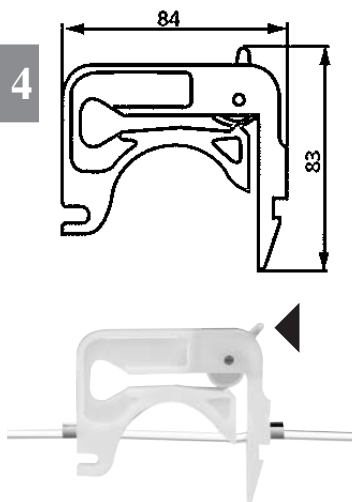
MS/CA Anpresshebel
MS/CA pressure lever
MS/CA levier de pression

3 Color Code Stopper Schläuche
3-stop colour-coded tubing
Tubes color-codé à 3 arrêts

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Type of cassettes

MS/CA Click'n'go

Spring-loaded cassette with stainless steel pressure spring providing defined and reproducible occlusion conditions. Independent of tubing diameter and material, the occlusion is automatically correct.

MS/CA Pressure lever

Cassette with pressure lever for adjustable occlusion (may be an advantage for elevated differential pressure conditions).

Cassette material

POM-C

(Polyoxymethylene-Copolymer) (Polyacetal, Polyformaldehyde) Good chemical resistance to many organic solvents and strong alkaline chemicals. Is affected by strong acids and oxidizing substances. UV-stabilized and stable at (dry) temperatures up to 80°C (in continuous use) or 136°C (for a short time).

PVDF ¹⁾

(Polyvinylidene fluoride) Very good chemical resistance to acids and most aliphatic, aromatic and chlorinated solvents. Not suitable for long contact with esters, ketones, amines and strong alkaline chemicals. Stable to UV and temperatures up to 110°C (continuous use) or 142°C (for a short time)

¹⁾ only cassettes with pressure lever

Pump tubing

These cassette pumps take tubing with 3 colour-coded collars. Thanks to these collars the tubes are always inserted under the same conditions, which results in a high repeatability of the flow rate. What is more, the different colour codes help to easily identify the various tubing diameters.

Tubing with 3 collars can be used twice. Thanks to this advantage the service life of the tubes can be doubled.

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